

IHC PROTOCOL

IHC standard protocol optimized for Triple A Polyclonals and Precisa Monoclonals from Atlas Antibodies.

DEPARAFFINIZATION

Paraffin sections of 4 µm thickness are baked overnight at 50°C. Prior to immunostaining, deparaffinization and hydration is performed in xylene and graded ethanol to distilled water.

During hydration, a 5 minutes blocking for endogenous peroxidase is performed in 0.3% H₂O₂ in 95% ethanol.

ANTIGEN RETRIEVAL

1. Standard antigen retrieval method

The standard antigen retrieval method is Heat Induced Epitope Retrieval (HIER) in retrieval buffer pH 6.1 (Dako, Agilent, Santa Clara, CA, USA) using a pressure boiler (Decloaking chamber NxGen, Biocare Medical, Pacheco, CA, USA) as heat source.

HIER is performed by heating the TMA-slides immersed in retrieval buffer for 20 minutes at 110°C in the pressure boiler. After completing the boiling, slides remain in the pressure boiler and are allowed to cool to 90°C. The total processing time is approximately 60 minutes

NOTE: The specified working dilutions of the primary antibodies are to be considered as a guideline only. Optimal dilutions must be determined by the user.

2. Alternative antigen retrieval method

For selected antibodies, alternative retrieval buffers and/or enzymatic antigen retrieval may be used as stated in the Product Datasheet and on the Antibody/ Antigen information page on the Human Protein Atlas website.

Enzymatic antigen retrieval method

Enzymatic retrieval is performed in the immunostaining instrument by incubating TMA-slides in Proteinase K solution (Lab Vision, Fremont, CA, USA) for 10 minutes at room temperature.

Heat-Induced Epitope Retrieval (HIER) in Retrieval Buffer pH 9

HIER is performed in Target Retrieval solution pH 9 (Dako, Agilent, Santa Clara, CA, USA) using Decloaking Chamber NxGen as described above.

IMMUNOHISTOCHEMICAL STAINING PROGRAM

All incubations are performed at room temperature.

All reagents are applied at a volume of 300 µl per slide.

1. Rinse in wash buffer.*
2. Incubate with Ultra V Block for 5 minutes.
3. Rinse in wash buffer (x2).
4. Incubate with primary antibody for 30 minutes.
5. Rinse in wash buffer (x3).
6. Incubate with Primary Antibody Enhancer for 20 minutes.**
7. Rinse in wash buffer (x2).
8. Incubate with labeled polymer for 30 minutes.
9. Rinse in wash buffer (x2).
10. Develop in DAB solution for 5 minutes.
11. Rinse in distilled water.
12. Counterstain in hematoxylin for 5 minutes.***
13. Rinse in tap water for 5 minutes.
14. Rinse in lithium carbonate (Li₂CO₃) water, diluted 1:5 from saturated solution for 1 minute.
15. Rinse in tap water for 5 minutes.
16. Dehydrate in graded ethanol and xylene
17. Mount in Pertex.
18. Coverslip.

* Steps 1 – 11 are performed in Autostainer 480S (ThermoFisher Scientific, Waltham, MA, USA).

** For polyclonal antibodies skip steps 6 and 7.

*** Steps 12 -18 are performed in fully automated integrated stainer Leica ST5010-CV5030 (Leica Biosystems Nussloch GmbH, Nussloch, Germany).

Warranty: The products supplied by Atlas Antibodies are warranted to meet stated product specifications and to conform to label descriptions when used and stored properly. Unless otherwise stated, this warranty is limited to one year from date of sales for products used, handled and stored according to Atlas Antibodies AB's instructions. Atlas Antibodies AB's sole liability is limited to replacement of the product or refund of the purchase price. All products are supplied for research use only. They are not intended for medicinal, diagnostic or therapeutic use. No products from Atlas Antibodies may be resold, modified for resale or used to manufacture commercial products without prior written approval from Atlas Antibodies AB Rev. December 2012

IHC PROTOCOL

REAGENTS

For immunohistochemistry, the following reagents are commercially available from Thermo Fisher Scientific, Waltham, MA, USA:

1. Wash buffer (10x concentrate). Working solution originally contains 0.05% (v/v) Tween 20. Extra Tween 20 is added to a final concentration of 0.20%.
2. Antibody diluent
3. UltraVision LP HRP polymer®
4. Primary Antibody Enhancer (only for monoclonal antibodies)
5. Ultra V Block
6. DAB Quanto Detection System (including chromogen and substrate)

For epitope retrieval, the following buffers are commercially available from Dako, Agilent, Santa Clara, CA, USA:

- Target Retrieval Solution, Citrate pH 6.1 (10x)
- Target Retrieval Solution, pH 9 (10x)

In addition, Mayer's hematoxylin and xylene (Histolab, Gothenburg, Sweden) are used.

Revision C, November 2021

Warranty: The products supplied by Atlas Antibodies are warranted to meet stated product specifications and to conform to label descriptions when used and stored properly. Unless otherwise stated, this warranty is limited to one year from date of sales for products used, handled and stored according to Atlas Antibodies AB's instructions. Atlas Antibodies AB's sole liability is limited to replacement of the product or refund of the purchase price. All products are supplied for research use only. They are not intended for medicinal, diagnostic or therapeutic use. No products from Atlas Antibodies may be resold, modified for resale or used to manufacture commercial products without prior written approval from Atlas Antibodies AB Rev. December 2012